



Performance Qualification

• Precision of the Stand. Gradient

• Instruments and Fluidics

Instrument Name	Model	Supplier's Name	Serial Number
Pump	LPG-3400A	Thermo Scientific	8002568
UV Detector	DAD-3000RS	Thermo Scientific	8043573
Chromeleon Datasystem	V. 6.80 SR16 Build 5387	Thermo Scientific	40640

Accessories	Description
Back Pressure Device	Capillary (L:15 m; ID:0,18 mm)
Solvent A	Water (HPLC-Grade)
Solvent B for Gradient	Water + 0.1% Acetone

• Additional Information

Customer: Customer's Name
 Operator: Operator's Name
 Operator's Jobtitle

Execution Date: 2020-mai-27
 Next Qualification: 2020-nov


• Limits and Test Results

	Limit	Observed max. SD	Result of all Steps
Gradient Precision	<= 0,500 %	0,080 %	Test passed

• Signatures

	Signature	Date
Submitter / Operator:	{seq.submitOperator.userName}	{seq.submitTime}
Reviewer:	{seq.reviewOperator.userName}	{seq.reviewTime}
Approver (e-sig. only):	{seq.approveOperator.userName}	{seq.approveTime}

• Overlay of three Gradients

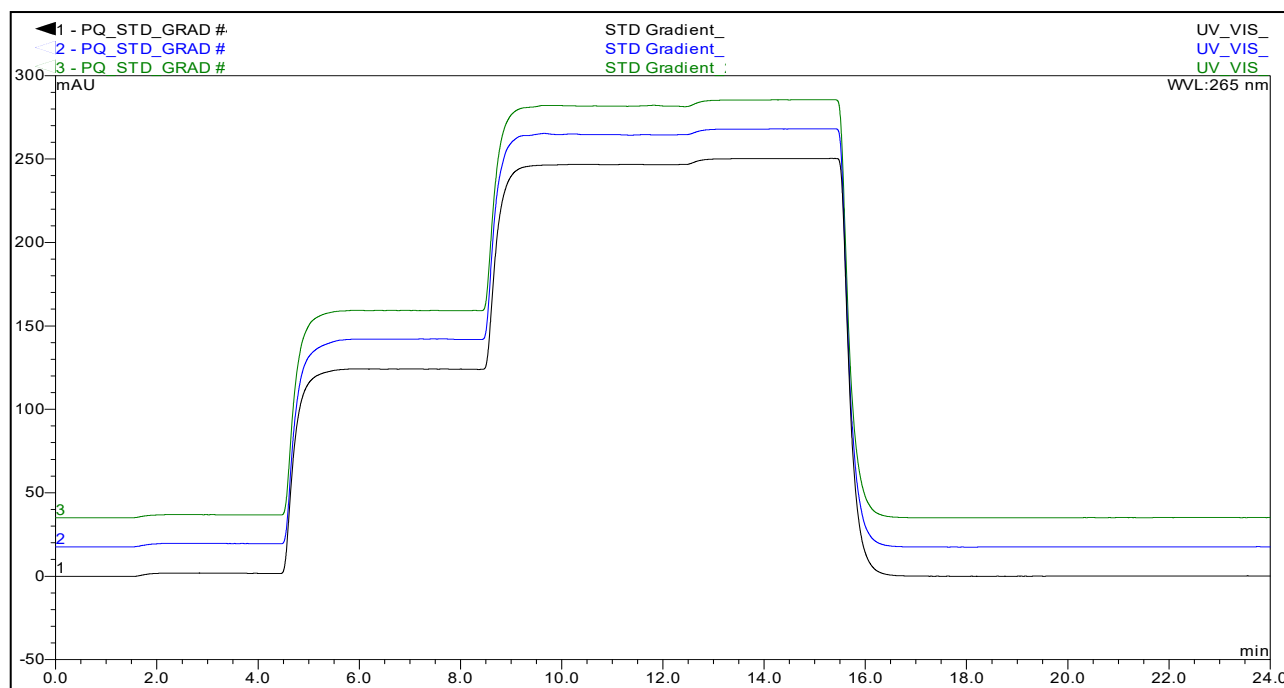

• Data of three Gradients

Expected Value [%]	Calculated Value Gradient_1 [%]	Calculated Value Gradient_2 [%]	Calculated Value Gradient_3 [%]	Critical SD [%]	Calculated SD [%]	Result
0,00	0,00	0,00	0,00	0,50	0,000	ok
1,00	0,81	0,70	0,70	0,50	0,063	ok
50,00	49,71	49,56	49,59	0,50	0,080	ok
99,00	98,59	98,64	98,56	0,50	0,040	ok
100,00	100,00	100,00	100,00	0,50	0,000	ok

Determination of Pump Unit for Dionex DGPs

Sequence name: PQ_STD_GRAD

Right end of the sequence GRAD

Chromeleon (c) 2017 Thermo Fisher Scientific Inc.
Version 6.80 SR16 Build 5387 (291597)

Gradient STD / PUMP_GRADIENT_REPRO
Printed: 29/05/2020 16:14

Pump's model number: LPG-3400A
Pump's model variant: LPG
Pump unit:

Determination Page Title

Gradient type:	• Precision of the Stand. Gradient
Solvent C-D (if used)	
Pump unit (DGPs only)	
All:	• Precision of the Stand. Gradient



